

The LabGard® ES (Energy Saver) AIR

# NU-543

# NUAIRE



environmental,  
product & personnel  
protection

The LabGard® ES (Energy Saver) AIR model NU-543 is a Class II, Type A2 Biological Safety Cabinet offering Personnel, Product, and Environmental Protection through the use of quality design, an ultra-high-efficiency motor, and longer lasting filters all at an affordable price.

You deserve a quality and robust Biological Safety Cabinet every time you sit down to work. Every NuAire Biological Safety Cabinet goes through a series of quality assurance checks before it reaches your laboratory. As your product moves through the assembly plant, each work station double checks the work of the previous station,

filters are triple-checked, and biological tests are continually being performed under dynamic conditions to ensure maximum safety.

Each cabinet is built to order for your application. It is as simple as including an additional service valve or as complex as developing the interior to fit a specific piece of machinery.

Our local sales representatives are the best in the industry. They assist with product selection, built-in options, and accessories, discuss pricing, offer installation guidance, and provide support over the life of your NuAire product.

# The LabGard® ES (Energy Saver) AIR NU-543




## Flexibility and Innovative Design

### Maximum Access Flexibility

Select from three window access opening options to meet the concerns of your facility and/or staff.

- **8"** (203 mm) lower operating cost (NSF/ANSI 49)
- **10"** (254 mm) lower costs and improved ergonomics (NSF/ANSI 49, EN12469)
- **12"** (305 mm) improved ergonomic reach by 40%\* (NSF/ANSI 49)

\* As compared to an 8" (203 mm) access opening.

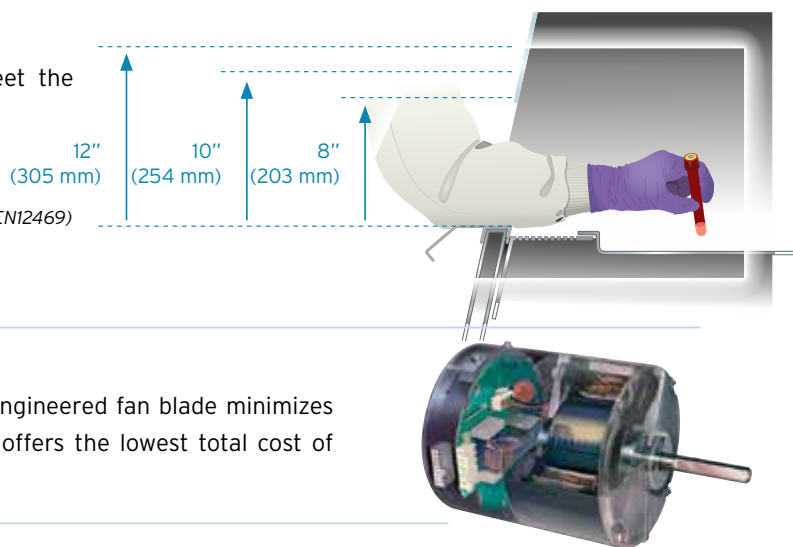

### Ultra High Efficiency Motor

A single ultra-high efficiency ECM motor with a precision engineered fan blade minimizes energy usage while extending the life of filters. LabGard® offers the lowest total cost of ownership compared to other models.

### A Size to Fit Your Facility or Application

3 FT (0.9 M)



4 FT (1.2 M)



5 FT (1.5 M)



6 FT (1.8 M)



# The LabGard® ES (Energy Saver) AIR NU-543




## Ergonomic Design

NuAire designs equipment for you that provides improved safety conditions; improved productivity; improved quality and reliability; and reduction of arm, shoulder, and neck strain which can contribute to common work place injuries.



- ADA Compliant
- More Usable Work Area
- Shorter Reach Into Work Zone
- Recessed Air Foil
- Recessed Work Zone
- More Knee Room
- Ability to Sit or Stand at a Range of Heights
- 21" (533 mm) Frame-less Edge Window Sash Improves Vision/Sight Lines Into Work Zone
- Staggered Service Valves
- Accessories: Elbow rests, Armrest Foam Pads, Ergo chair, Adjustable Base Stand, Stainless Steel Turntable

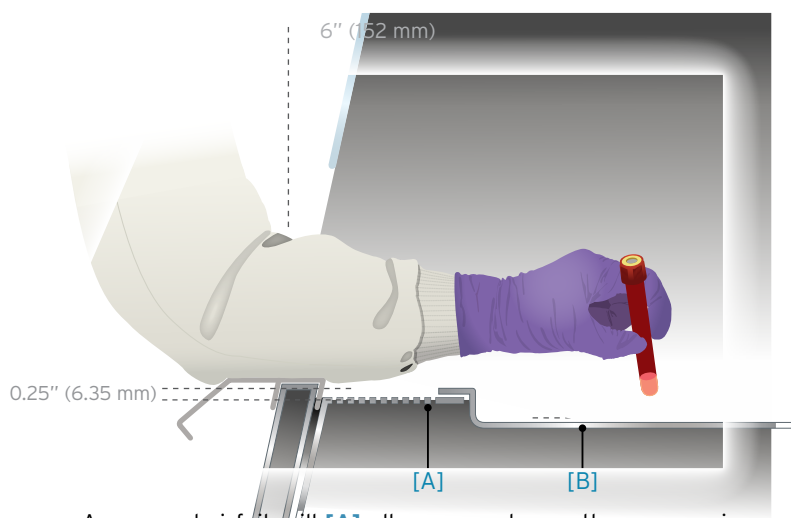


A recessed airfoil grill [A] allows users to use the ergonomic armrest without blocking airflow compromising safety. A recessed work tray [B] provides users greater flexibility when working with bottles or pipettes. A shorter reach into the work zone allows users to bring work closer without sacrificing safety. These are just some of the user focused ergonomic designs NuAire® incorporates into all our laboratory products.

The LabGard® ES (Energy Saver) AIR

# NU-543

**NUAIRE**

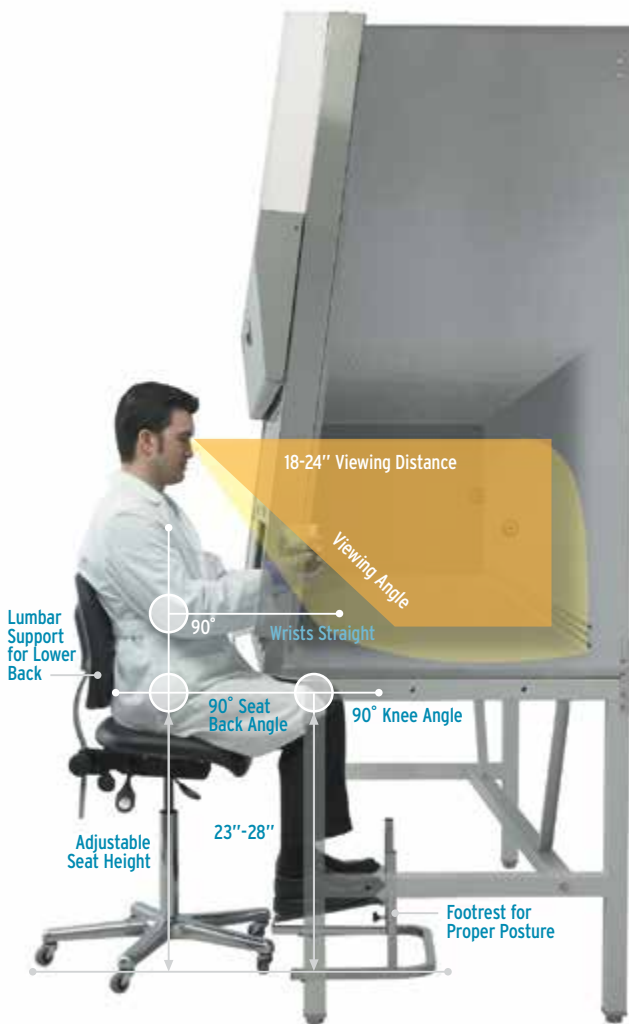

[ Larger Work Area ]



[ Expanded Vision Zone ]



[ Disposable Arm Rest Pads ]



[ Recessed Airfoil Grill and elbow rest accessory ]

# The LabGard® ES (Energy Saver) AIR NU-543




## Quality - Nothing is Built like a NuAire

**Monolithic Design.** At core of any biosafety cabinet is its basic structure. NuAire has over 100,000 biosafety cabinets in the field; some of which have been running strong since the 1970's.

The NuAire design is a pressure tight, single piece of type 304 stainless steel. A stainless steel wrap [A] makes up the walls of the cabinet. Stainless steel sheets [B] are seam welded in place to make up the top and bottom of the cabinet. A smaller wrap [C] featuring coved interior corners is seam welded in place to make up the back and side walls of the cabinet. This improves the cabinet integrity by creating a double sidewall. Work tray supports [D] help minimize the vibration transfer from the shell of the cabinet to the work zone.


The end result is a pressure tight monolithic shell of 100% stainless steel [E] without the use of silicone like many multi-piece cabinets. This provides you with a solid foundation for longer product life eliminating potential leaks.

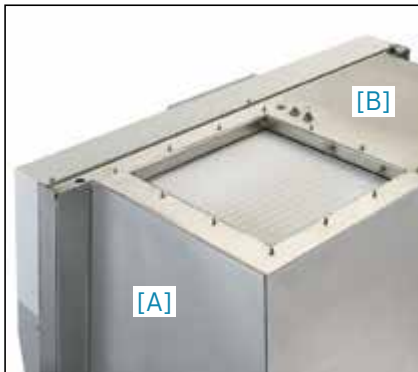

The LabGard® ES (Energy Saver) AIR

# NU-543

**NUAIRE**

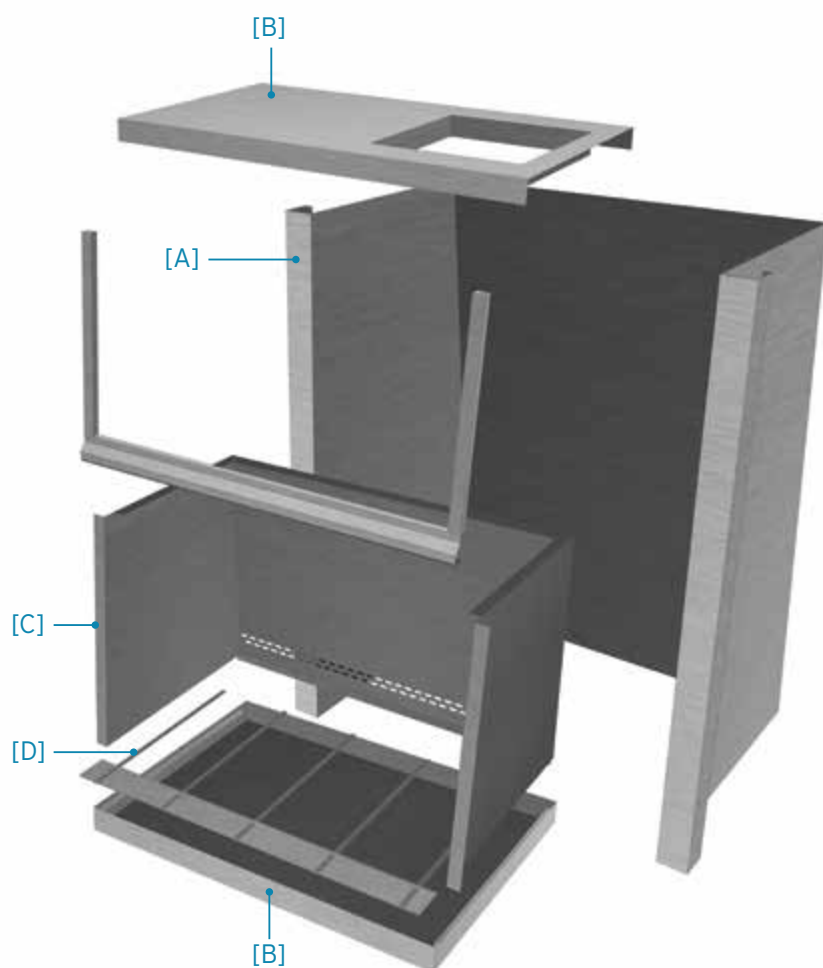




## Efficient, Quiet, Powerful

The LabGard® AIR comes standard with an Ultra High Efficiency Motor. Each motor has a precision engineered blower fan to maximize energy efficiency and filter life. The more powerful motor in model NU-543 equates to longer filter life lowering your total cost of ownership.

### Power Savings

Model	Nominal Width	Motor Power	Savings*
NU-543-300	3 ft (0.9 m)	¾ HP	59%
NU-543-400	4 ft (1.2 m)	¾ HP	
NU-543-500	5 ft (1.5 m)	1 HP	
NU-543-600	6 ft (1.8 m)	1 HP	

\* Compares NuAire DC ECM motor technology and a BSC featuring DC motor technology. Information taken from page 7.

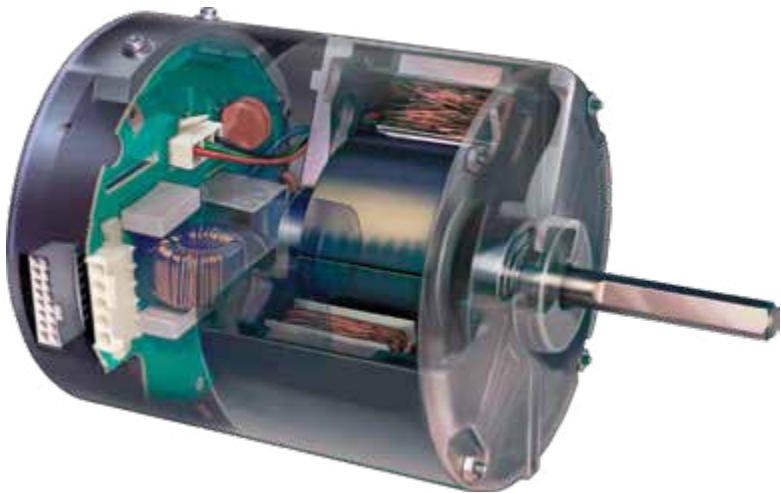

## Your Safety is Our Number One Priority

NuAire's commitment to safety doesn't stop at meeting industry standards. NuAire continually performs dynamic tests to push the limits of our product's performance to ensure you and your laboratory personnel will remain safe.

Automatic Airflow Compensation - the Ultra High Efficiency motor will automatically increase or decrease speed to maintain proper airflow settings based on readings from FlowGard Electronic Control System.



True laminar (unidirectional) downward moving air flushes the work surface at 60 fpm (0.30 m/s) minimizing cross contamination.

**true**laminar



A dynamic air barrier at the front of the cabinet separates the laboratory from the work zone environment at an air speed of 105 fpm (0.53 m/s).

*dynamic*  
**air**  **barrier**



The HEPEX™ Zero Leak Airflow system evenly disperses air over the supply and exhaust HEPA filters, minimizes vibration transfer, and eliminates potential leaks by surrounding positive pressure chambers with negative pressure.

**HEPEX**  **zero-leak**  
**airflow**




## Make it Easier on your Service Personnel

We at NuAire® have designed each LabGard® Biosafety Cabinet to be service friendly. **100%** of service functions can be performed at the front or top of the cabinet.

### Drop Down Control Panel

A flip down control panel creates a workable tray for service technicians.

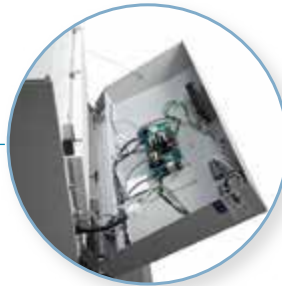

### Exhaust Choke

Easy access to exhaust choke to certify airflows.




### Front Filter Access

Access filters and the motor blower from the front of the cabinet with the removal of a safety panel.




### DECON101 System

Easily connect to a variety of decontamination systems that use Formaldehyde, Chlorine Dioxide or Vapor Phase Hydrogen Peroxide

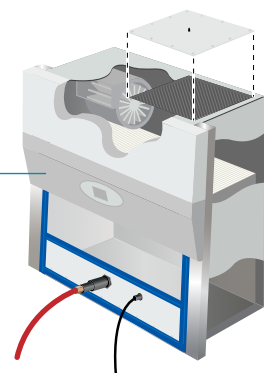

Controlled Safety - **FLOWGARD** Electronic Control System

The LabGard® AIR offers the latest digital microprocessor design technology for improved cabinet performance and safety. The FlowGard™ control system integrates a dual thermistor airflow sensor (*intelliflow*™) to monitor the cabinet's airflow performance. The FlowGard™ control system also integrates an ultra high efficiency motor / blower that provides automatic compensation for both filter loading and line voltage variances. The FlowGard™ also controls the motor / blower, visible light, UV light (optional), outlet(s), alarms, and the sliding window position via micro switches.

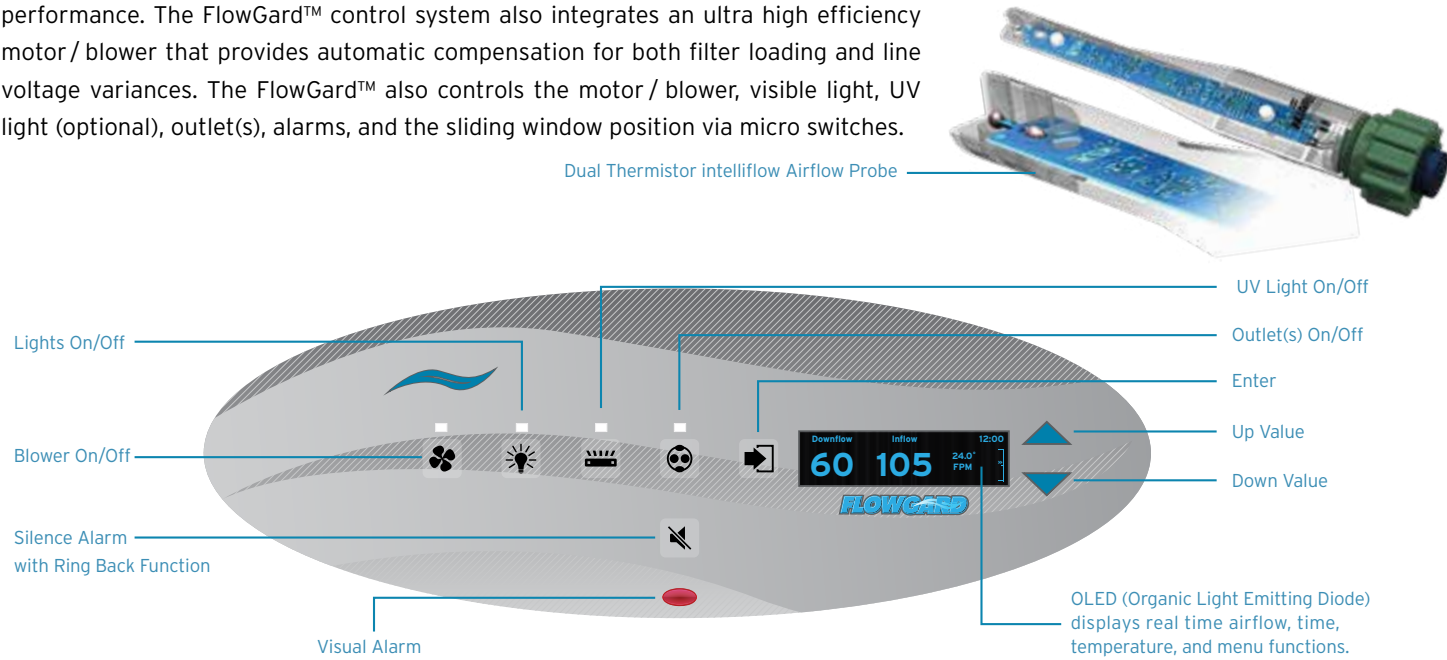


**nitecare**

Night Setback Mode - Upon sliding the window sash into the closed position, the motor blower will continue to operate at a lower rate to save energy and maintain interior clean air conditions ready for use upon window sash opening.

## Perfectly Designed for You and Your Safety

### Airflow Protects Personnel and Product

LabGard® AIR utilizes a single motor blower system to maintain airflow performance. The single motor blower system is inherently balanced. As particulates load the HEPA filters between certifications, NuAire's single motor/blower system maintains a consistent ratio of down flow to inflow, preventing uneven filter loading. This maintains a safer environment for you and your laboratory as well as maintains the proper down flow to minimize cross contamination within the work zone.

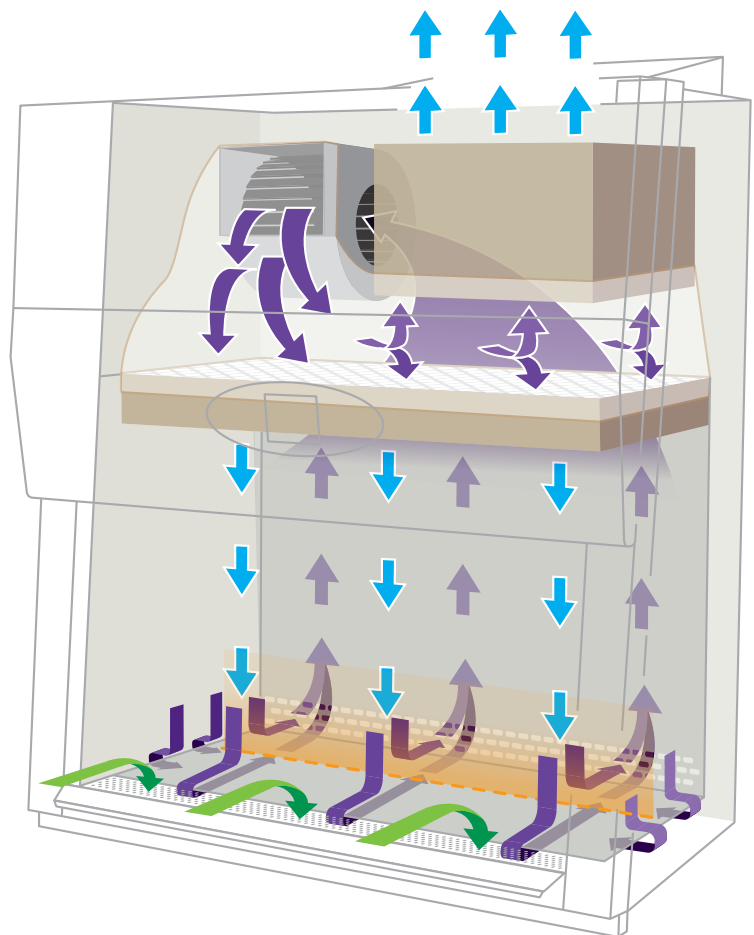

# The LabGard® ES (Energy Saver) AIR NU-543






The LabGard NU-543 AIR is a Class II Type A2 Biological Safety Cabinet that can be either exhausted back into the room or connected to a facility HVAC system. The cabinet's airflow is 30% exhausted / 70% recirculated to minimize cross-contamination of sterile hazardous drugs and low to moderate risk biologicals in the absence of volatile toxic chemicals.

## Specifications

Model	Width Size (nominal)	Electrical (all sizes)	Interior Dimensions (W x D x H) in [mm]	Exterior Dimensions (W x D x H) in [mm]	Net Weight
NU-543-300	3 ft [ 0.9 m ]	115V 60Hz E/S: 230V 50/60Hz K: 100V 60HZ J: 100V 50/60Hz	34 <sup>3</sup> / <sub>8</sub> x 25 <sup>3</sup> / <sub>4</sub> x 28 <sup>1</sup> / <sub>2</sub> [ 873 x 654 x 724 ]	41 <sup>5</sup> / <sub>8</sub> x 31 <sup>7</sup> / <sub>16</sub> x 60 <sup>7</sup> / <sub>8</sub> [ 1057 x 799 x 1546 ]	410 lbs [ 186 kg ]
NU-543-400	4 ft [ 1.2 m ]		46 <sup>3</sup> / <sub>8</sub> x 25 <sup>3</sup> / <sub>4</sub> x 28 <sup>1</sup> / <sub>2</sub> [ 1178 x 654 x 724 ]	53 <sup>5</sup> / <sub>8</sub> x 31 <sup>7</sup> / <sub>16</sub> x 60 <sup>7</sup> / <sub>8</sub> [ 1362 x 799 x 1546 ]	480 lbs [ 218 kg ]
NU-543-500	5 ft [ 1.5 m ]		58 <sup>3</sup> / <sub>8</sub> x 25 <sup>3</sup> / <sub>4</sub> x 28 <sup>1</sup> / <sub>2</sub> [ 1483 x 654 x 724 ]	65 <sup>5</sup> / <sub>8</sub> x 31 <sup>7</sup> / <sub>16</sub> x 60 <sup>7</sup> / <sub>8</sub> [ 1669 x 799 x 1546 ]	570 lbs [ 258 kg ]
NU-543-600	6 ft [ 1.8 m ]		70 <sup>3</sup> / <sub>8</sub> x 25 <sup>3</sup> / <sub>4</sub> x 28 <sup>1</sup> / <sub>2</sub> [ 1788 x 654 x 724 ]	77 <sup>5</sup> / <sub>8</sub> x 31 <sup>7</sup> / <sub>16</sub> x 60 <sup>7</sup> / <sub>8</sub> [ 1972 x 799 x 1546 ]	640 lbs [ 290 kg ]

\* Specify models with appropriate letter suffix for electrical specifications. "NU-543-400E" for 230 VAC 50/60 Hz

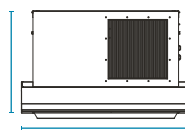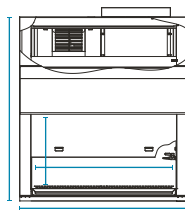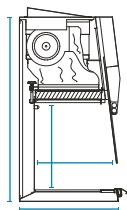

The LabGard® ES (Energy Saver) AIR

NU-543



Features

Performance

Airflow 30% exhausted / 70% recirculated  
Dynamic Air Barrier 105 fpm (0.53 m/s) Inflow  
True Laminar Airflow 60 fpm (0.30 m/s) Down Flow  
Filter Load Capacity 300% Plus  
Filter Life ±10 Years  
NSF/ANSI 49, EN12469

Protection

HEPEX Zero Leak Airflow System  
ANSI/ASHRAE 110, Surrogate Powder, Biological Tested and Validated  
Airflow / Window Alarms with Ring Back Function  
Security Password Protection  
UL Listed, CE

Efficiency

Ultra High Efficiency ECM Eon™ Motor with Precision Engineered Forward Curved Fan  
Auto Filter Loading Compensation

Motor Speed Controller  
Internal Exhaust Damper  
nitecare Night Setback

Construction

Monolithic 16/18 Gauge, Type 304 Stainless Steel  
Welded Pressure Tight Construction (Silicone Free)  
Coved Interior Corners  
Removable Control Center  
Removable Stainless Steel Coved Work Tray  
Recessed and Removable Front Intake Grill  
Stainless Steel Spill Trough with Drain Valve  
Vibration Reducing Work Surface Support System  
External Non-Glare 100 lm/ft<sup>2</sup> Fluorescent Lighting  
Frame-less Sliding Viewing Window  
Metal Diffuser Over Supply Filter  
Permanent Positive Pressure Plenum with Quick Release Supply Filter Removal

Easy transport design  
100% Front Service Friendly

Ergonomics

Comfort 10° sloped front window  
ADA Compliant  
Large User-friendly Work zone with coved corners (20-3/4 inch (527mm) depth)  
Low noise ±62 dba operation with Ambient Conditions of 53 dba  
Maximum visibility into work zone 23.75-inches (603 mm) viewing area  
Easy slide front window sash, 21-inches (533 mm) full open  
Centrally located control panel  
Attenuomont Vibration Control

Design Your Own

Nominal Width

3-foot (0.9 m)  
4-foot (1.2 m)  
5-foot (1.6 m)  
6-foot (1.8 m)

Access Opening

12" [305 mm] Access Opening (NSF/ANSI 49)  
10" [254 mm] Access Opening (NSF/ANSI 49, EN12469)  
8" [203 mm] Access Opening (NSF/ANSI 49)

Supply Filter

HEPA Filter 99.99% efficient @ 0.3 μ (NSF/ANSI 49)  
HEPA H14 Filter 99.995% efficient @ 0.025 MPPS (EN12469)

Exhaust Filter

HEPA Filter 99.99% efficient @ 0.3 μ (NSF/ANSI 49)  
HEPA H14 Filter 99.995% efficient @ 0.025 MPPS (EN12469)  
Redundant (Qty 2) HEPA H14 Filter 99.995% efficient @ 0.025 MPPS (EN12469)

Power

Power Cord  
Junction Box

External Lighting

Fluorescent [90 - 120 fc (968 - 1291 lux)]  
LED [90 - 120 fc (968 - 1291 lux)]

Armrest (Ergonomic)

Heavy duty plastic  
Metal

Options

Ultraviolet Light  
IV Bar with 3 Height Locations, 6 Hooks  
NSF Approved Cord Pass Through  
Stainless Steel Window Edge Protector

Prop-up Work Tray  
Australian Test Method

Sidewalls

Stainless Steel  
Glass

Service Valves/Couplings

Smooth Interior  
Left/Right Bottom  
Left/Right Middle  
Left/Right Top

Plumbing

Soft Copper Pipe 3/8"  
Black Pipe 3/8" (Gas Only)  
Rigid Copper 1/2"

Decorative Side Panels

None  
No Plumbing  
Plumbing Out Top  
Plumbing Out Back  
Plumbing Out Bottom

Outlets (Quantity 2)

Duplex Outlet (NEMA 5-15) (3 Amp Max)  
GFI Duplex Outlet (NEMA 5-15) (3 Amp Max)  
GFI Duplex Outlet (NEMA 5-15) (5 Amp Max)  
Duplex Outlet (NEMA 5-15) (15 Amp Max)  
GFI Duplex Outlet (NEMA 5-15) (15 Amp Max)  
Duplex Outlet (NEMA 5-20) (20 Amp Max)  
GFI Duplex Outlet (NEMA 5-20) (20 Amp Max)  
Duplex Outlet (NEMA 6-15) (15 Amp Max)  
Type D - United Kingdom  
Type E - France  
Type F - Schuko  
Type G - United Kingdom

Type H - Israel  
Type I - Australia  
Type J - Switzerland  
Type K - Netherlands  
Type M - South Africa

Seismic Brackets

Side Mount  
Rear Mount  
Floor Mount

Custom Solutions

Microscope Window  
Sinks with Hot/Cold or DI water Faucets  
Heated Work-surface  
Widths and/or Heights

Base Stands

Telescoping Base Stand with Leg Levers or Casters  
Automatic Adjustable Base Stand with Leg Levers or Casters  
Base Storage Telescoping Base Stand with Leg Levers or Casters  
Base Stand Shelf [8 or 12-inch width (203 or 305 mm)]  
Ergo-tron Workstation for Monitor and Keyboard

Exhausting

Variable Range Low Flow Canopy (VFC) with Monitor  
External Airflow Alarms  
Flex Duct Kits  
Butterfly Dampers  
Silicone Connection Sleeve Kit  
Round to Round Transition [8 to 10-inch (203 to 254 mm)]

Ergonomics

Adjustable Footrest  
Stainless Steel Turn Table [12-inch (305 mm) Diameter]  
Small Elbow Rest [3 x 5-inches (76 x 127 mm)]  
Large Elbow Rest [3 x 10-inches (76 x 254 mm)]  
Foam Armrest Pads



# The LabGard® ES (Energy Saver) AIR NU-543



Each laboratory or pharmacy is unique and every application can differ. Set your laboratory up for success by designing your own NuAire Biosafety Cabinet. It can be as simple as adding a service valve or as involved as customizing the height of the work zone. NuAire can do it all. Invest in a product that lasts and is designed for your needs.

## Popular Configurations:

### Configuration 1 - Hazardous Drug Compounding



Smooth Interior for clean-ability



IV Bar with Three Locations and Six Hooks



Cord Pass-through (NSF Approved)



Stainless Steel Armrest



Prop Up Work Surface



VFC Exhaust Transition with Alarm



Silicone Sleeve Exhaust Connection Kit



Telescoping Base Stand with Ergotron Computer Monitor and Keyboard Arm and Base Stand Shelf

### Configuration 2 - Tissue Culture



Service Valves/Couplings [Gas, Air, Vac]



Ultraviolet (UV) Light



Cord Pass-through (NSF Approved)



Prop Up Work Surface



Motorized Base Stand - Leg Levelers or Casters

The LabGard® ES (Energy Saver) AIR

# NU-543



### Base Stands



Telescoping Base Stand  
- Leg Levelers or Casters



Motorized Base Stand  
- Leg Levelers or Casters



Storage Telescoping Base Stand  
- Leg Levelers or Casters



Base Stand Shelf  
- 8 or 12-inch width (203 or 305 mm)



Ergotron® Arm for Monitor and Keyboard / Mouse

### Exhausting



VFC Exhaust Transition



Fixed Exhaust Canopy Transition w/ Alarm



Silicone Sleeve



Flex Duct



Manual and Automatic Butterfly Valves

### Ergonomics



Stainless Steel Turntable  
- [12-inch (305 mm) Diameter]



Elbow Rest  
- Small Elbow Rest [3 x 5-inches (76 x 127 mm)]  
- Large Elbow Rest [3 x 10-inches (76 x 254 mm)]




Adjustable Foot Rest

### Built-in Options



LED Lighting (Replaces Florescent Lighting)

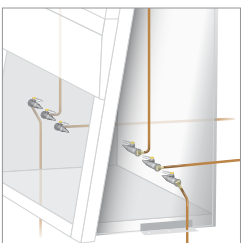


Pre-Piped Plumbing



Outlets



Cord Pass-through (NSF Approved)

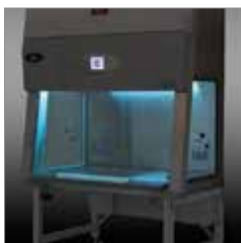

Ultraviolet (UV) Light



The LabGard® ES (Energy Saver) AIR  
**NU-543**



## Complete your Laboratory

NuAire offers a variety of Laboratory Equipment to meet the needs of the research, vivarium, pharmacy, environmental laboratory and more. Join the NuAire family and complete your laboratory with complementary NuAire Quality products.



**LABGARD**  
Biosafety Cabinets



**IN-VITROCELL<sup>ES</sup>**  
CO<sub>2</sub> Incubators



**NUWIND**  
Centrifuges



**Koki**  
CENTRIFUGES



**BLIZZARD**  
Ultralow Freezers



**AIRECARD<sup>ES</sup>**  
Laminar Airflow Workstations



**PHARMAGARD<sup>ES</sup>**  
Compounding Isolators



**ALLERGARD<sup>ES</sup>**  
Animal Laboratory Products



**Polypropylene Fume Hoods  
and Casework**



**Custom Solutions**

# The LabGard® ES (Energy Saver) AIR NU-543



**Delivering product to the customer that is cost effective, exceeds their requirements, delivers product on-time, and increases productivity and profitability.**

In every NuAire product you'll find brilliant but practical design, keen attention to detail in every phase of the fabrication and assembly process, thoroughly tested, outstanding value, and dependable service. These are the foundation for NuAire's international reputation, and practical reasons why NuAire is universally recognized as the finest. NuAire's Quality Policy to Satisfy Customers, Comply with Quality System Requirements and Continually Improve.

**Sales and Service Like No Other** NuAire employs a network of knowledgeable sales and service individuals some of which have been selling and servicing NuAire Laboratory Products since our inception in 1971.

**Working Safely Video Series.** Learn how to work safely in your NuAire Biological Safety Cabinet with this 7 Chapter online video series.



Our Sales Representatives are located throughout the globe providing you with a local individual who can assist product selection, pricing, installation, training, and service throughout the life of your product. Once you buy a NuAire you're a part of our family and we'll work hard to satisfy your needs.

NuAire offers service-training classes and individualize assistance to service professionals in the laboratory community. These NuAire certified companies are located throughout the world allowing NuAire to provide service and assistance to the farthest corners of the earth. Have the option to select from multiple service vendors in your area to find a company that fits your needs.

**Contracts, GPO's and many more!** NuAire has a variety of Contracts set up through Group Purchasing Organizations, Government Associations, Health Systems, Universities and many more.



ISO 9001:2015 Certificate

